



**Owner's Manual for the
PulseBlasterPlus! Spin Decoupler™**



**SpinCore Technologies, Inc.
3525 NW 67th Avenue
Gainesville, Florida 32653, USA
Phone: (352)-271-7383**

<http://www.spincore.com>

Congratulations and THANK YOU for choosing a design from SpinCore Technologies, Inc. We appreciate your business. At SpinCore we try to fully support the needs of our customers, so if you ever need assistance please contact us and we will strive to provide the necessary help.

© 2000-2002 SpinCore Technologies, Inc. All rights reserved. SpinCore Technologies, Inc. reserves the right to make changes to the product(s) or information herein without notice. PulseBlasterDDS™, PulseBlaster™, PulseBlasterPlus! Spin Decoupler™, SpinCore, and the SpinCore Technologies, Inc. logo are trademarks of SpinCore Technologies, Inc. All other trademarks are the property of their respective owners.

SpinCore Technologies, Inc. makes every effort to verify the correct operation of the equipment. This equipment should NOT, however, be used in system where the failure of a SpinCore device will cause serious damage to other equipment or harm to a person.

Contents

Section I: PulseBlasterPlus! Spin Decoupler™

1	Product Overview	3
2	System Architecture	4
3	User Guide	6
4	Connector Information	7

Appendix:

1	Available Options	8
---	-------------------	---

Section I: PulseBlasterPlus! Spin Decoupler™

1. Product Overview

The PulseBlasterPlus! Spin Decoupler™ system is an autonomous stand-alone pulse Radio-Frequency (RF) generator customized for NMR/MRI decoupling applications. The system integrates a precision timing engine and a dedicated Direct Digital Synthesis (DDS) core for generating continuous RF sequences for bi-level decoupling. To simplify the pulse-sequence programming, WALTZ and MLEV decoupling sequences are hard-coded via firmware and can be selected with an external TTL signal. The RF output frequency and pulse lengths can be controlled via a simple ASCII RS232 serial interface. The rf output power is approx. 10 dBm. The system is housed in a sturdy, shielded plastic enclosure with an integrated power supply.



Fig. 1 – A photograph of the PulseBlasterPlus! Spin Decoupler™

2. System Architecture

Figure 2 presents the general architecture of the PulseBlasterPlus! Spin Decoupler™ system. The four major sections of the design are as follows:

- Direct Digital Synthesis (DDS) core
- Precision Pulse Sequence Digital Timing (PP) core
- RF output (analog) circuitry, and
- RS-232 serial programming interface.

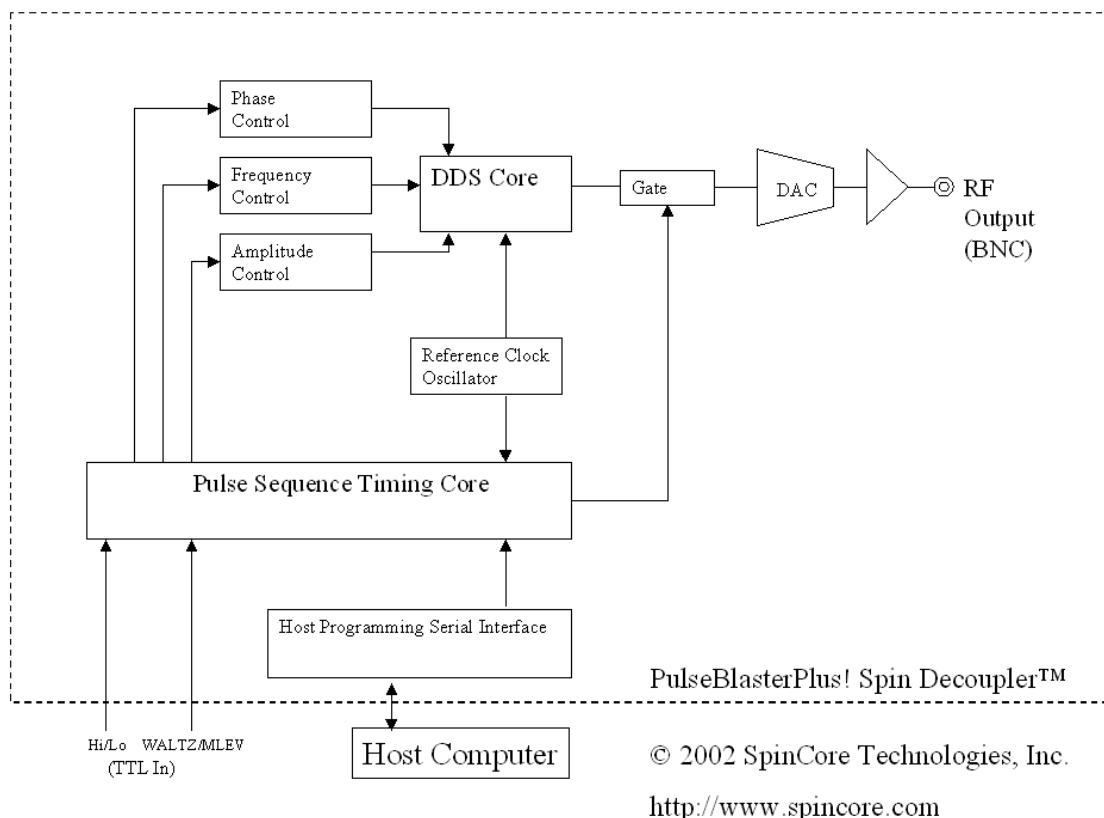


Fig. 2. PulseBlasterPlus! Spin Decoupler(tm) board architecture.

The DDS and PP cores run off a single, on-board reference clock oscillator. With a default 50 MHz clock oscillator, the DDS core allows for frequencies from 1kHz to 25MHz to be output, with a frequency resolution of 1kHz; the pulse-timing core allows for delay lengths of 1μs to 120ms, with a resolution of 1μs

The internal DDS core is a versatile digital frequency synthesizer with phase, frequency, and amplitude control registers to accommodate for broad requirements of RF pulse sequences used in NMR/MRI. A specific pulse decoupling sequence is produced via sequential control words generated

PulseBlasterPlus! Spin Decoupler™

by the pulse-sequence timing core. The timing core is an implementation of SpinCore's PulseBlaster™ pulse program timing processor - a dedicated, programmable microcontroller based on a 80-bit Very Long Instruction Word (VLIW) architecture. To simplify pulse-sequence programming, pulse programs are hard-coded in an EPROM. Two most-commonly used decoupling sequences, WALTZ and MLEV, have been programmed as default sequences. An external TTL signal (level) can select one of the two sequences. A second external TTL signal (level) can select between the full RF output power and a -3 dB value.

The RF output section converts the digital samples generated by the DDS core to analog waveforms, and it includes a high-speed Digital-to-Analog Converter (DAC). The gate between the DDS core and the DAC is continuously opened for the WALTZ/MLLEV decoupling sequences. The analog RF signals from the output of the DAC are amplified by a high-bandwidth amplifier delivering approx. 10 dBm of power; the output impedance is 50 ohm. The analog output signal is available on the external BNC connector.

The host programming serial interface, implemented as an RS-232 circuit operating at the fixed speed of 115 kBaud, allows the user to modify the output frequency and the pulse length of the generated pulse decoupling sequences. Any terminal/modem software capable of communicating via RS-232 communication channels (8-bit data, no parity, one stop bit) can be used to communicate with the system.

The system is NOT equipped with an interpolating filter. Thus, a step-like output signal will be observed, especially pronounced at high output frequencies. This is normal for non-filtered DDS-generated signals. To eliminate the step-like output, the user should use a low- or band-pass filter as appropriate for their application.

The system boots with the following parameters:

- Output frequency – 6.25 MHz
- Pulse sequence – WALTZ
- 90-degree pulse length – 2 ms
- RF output level – high (approx. 10 dBm).

By default, the system runs continuously once the board is powered.

3. User Guide

The PulseBlasterPlus! Spin Decoupler™ system is a stand-alone unit and is ready to use right out of the box. Upon receiving and unpacking the system, examine it for any damages that may have occurred in transit.

Attach a 50 ohm coaxial cable to the BNC output connector mounted on the box. For proper operation, the cable should be terminated with a 50 ohm impedance. Output signals can be evaluated with an oscilloscope with a 50 ohm input impedance. The observed signals will be in the range of approx. 3Vpp. Use an external filter, if desired, to eliminate the step-like output.

The PulseBlasterPlus! Spin Decoupler needs only to be powered. Connecting to a serial port on a host computer will allow the user to change the frequency and pulse length of the pulse sequence by a simple ASCII terminal interface.

To control the system using the serial interface, follow the steps below:

- 1) Configure your serial port or terminal software to use a 115200 bits per second transfer speed, 8 data bits, no parity bits, 1 stop bit, and no flow control.
- 2) Turn on the power switch of the PulseBlasterPlus! Spin Decoupler™.
- 3) Your terminal program should display the SpinCore Technologies' logo. At this point you are ready to begin controlling the frequency and pulse length of the sequence.
- 4) Press 'D' (*note this must be a capital letter) to change the value of the delay register. This value should be entered in μ s, should not contain any commas, and the valid range of values is 1 to 120000. If you enter too many digits or an invalid value, the delay register will not be updated correctly.
- 5) Press 'F' (*note this must be a capital letter) to change the value of the frequency register. This value should be entered in kHz, should not contain any commas, and the valid range of values is 1 to 25000. If you enter too many digits or an invalid value, the frequency register will not be updated correctly.
- 6) To control the amplitude of the output signal, a TTL signal should be connected to the RCA connector labeled "Hi/Lo." A high TTL level signifies full-swing output, and a low TTL level signifies half-swing output. This connector has an internal pull-up resistor, so the default mode of operation is full-swing output voltage.
- 7) To control the sequence being used, a TTL signal should be connected to the RCA connector labeled "WALTZ16/MLV16." A high TTL level signifies the WALTZ sequence, and a low TTL level signifies the MLEV sequence. This connector has an internal pull-up resistor, so the default mode of operation is WALTZ.

PulseBlasterPlus! Spin Decoupler™

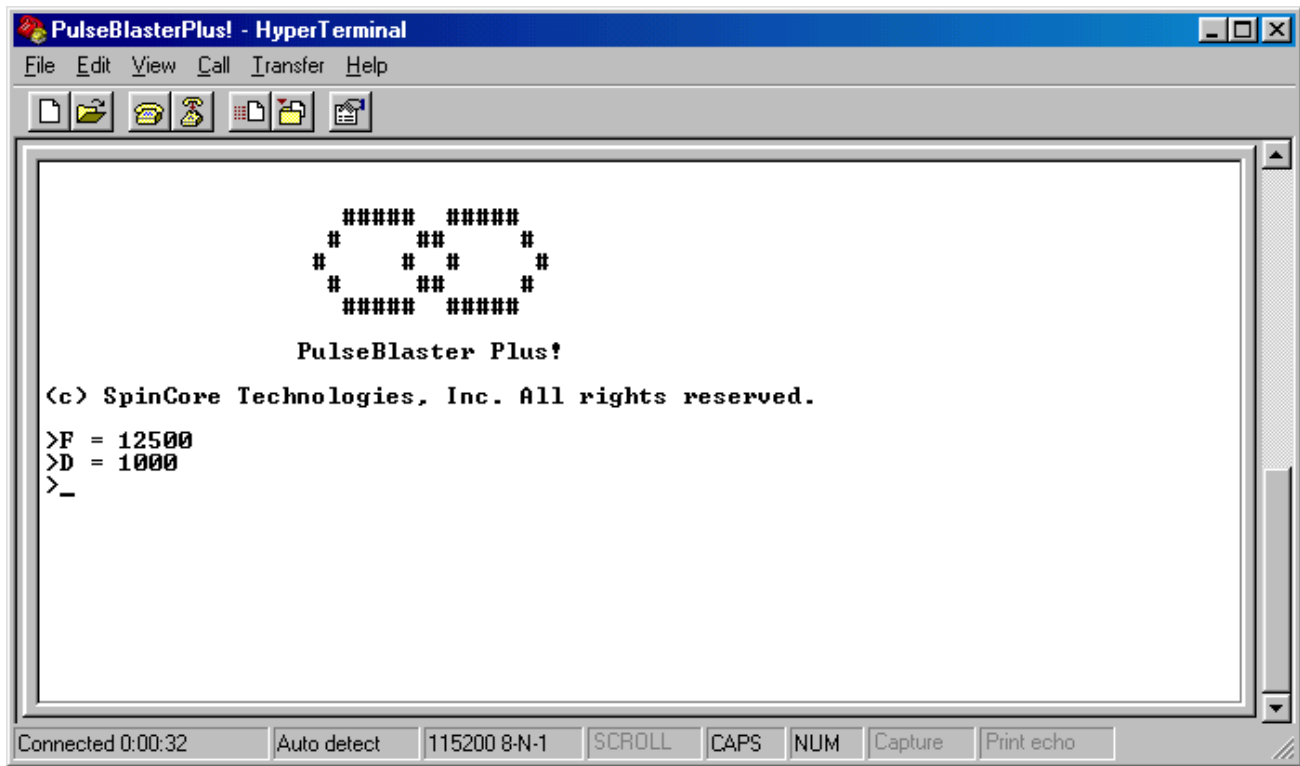


Fig. 3 – Microsoft Windows' HyperTerminal software can be used to control PulseBlasterPlus! Spin Decoupler™. Terminal settings: 115 kBaud, 8N1.

4. Connector Information

All input-output connectors are mounted on the back of the PulseBlasterPlus! Spin Decoupler™. A sketch depicting the available connectors is presented in Figure 4, below.

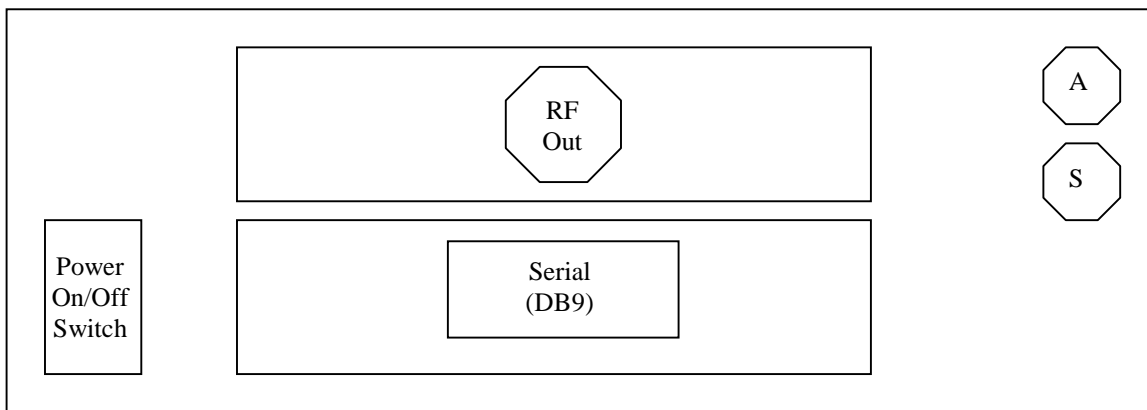


Figure 4 – Back view of PulseBlasterPlus! Spin Decoupler™

- RF Out – BNC connector for RF pulse sequence output signal; 50 ohm impedance, approx. 10 dBm.
- Serial (RS232 115kb) – DB9 connector to host computer serial port.
- A – RCA Jack Input TTL Level Amplitude Control (labeled Hi/Lo).
- S – RCA Jack Input TTL Level Sequence Control (labeled WALTZ/MLEV).

Appendix

1. Available Options

The PulseBlasterPlus! Spin Decoupler™ system can be customized. The following is a sample list of available options:

1. Reference clock frequency – any custom value of up to 100 MHz.
2. External gating.
3. Other pulse decoupling sequences.
4. Up to 8 additional TTL signals synchronized with the rf pulses.
5. Internal interpolating filter.
6. Internal power amplifier 30 dBm.
7. External power amplifier.

Please contact SpinCore Technologies, Inc., <http://www.spincore.com>, or sales@spincore.com, for more information.